


Microscope Manners

1. _____ carry a microscope _____ and with _____ hands- one hand on the _____ and one hand holding the _____.
2. Always return the _____ power objective to the position where it is pointing down to the _____.
3. Always bring the object into focus with the _____ power objective with the _____ knob and then use the _____ knob. Then move to the other _____ only using the _____ knob.
4. Never use the _____ focus knob to focus the _____ objective.
5. Never _____ the _____.
6. Don't _____ the microscope across the _____. _____ it up!

Microscope Manners

1. Always carry a microscope carefully and with 2 hands- one hand on the arm and one hand holding the base.
2. Always return the low power objective to the position where it is pointing down to the stage.
3. Always bring the object into focus with the low power objective with the coarse focus knob and then use the fine focus knob. Then move to the other objectives only using the fine focus knob.
4. Never use the coarse focus knob to focus the high power objective.
5. Never remove the eyepiece.
6. Don't drag the microscope across the counter. Pick it up!

Microscope Parts

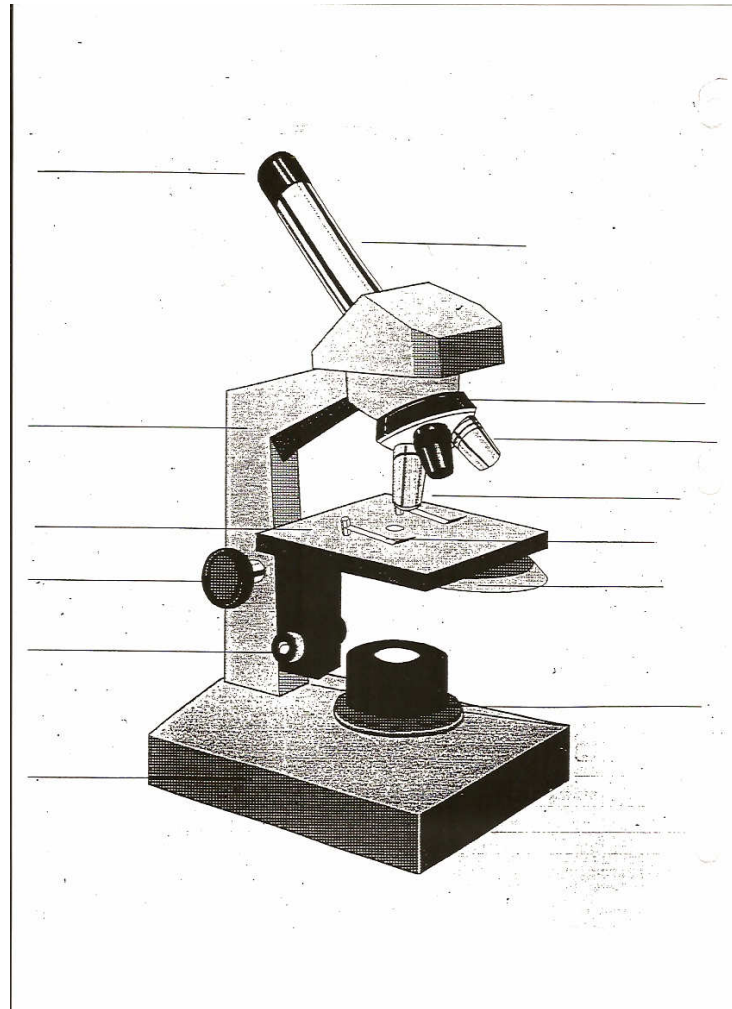

Down the left:
Eyepiece
Arm
Stage
Coarse Focus Knob
Fine Focus Knob
Base

Down the right:
Neck
Nosepiece
Low Power Objective
Stage Clips
Diaphragm
Light